

GingisKHAN®

FOR RESEARCH
USE ONLY
www.genovis.com

STORE AT
-20°C



SmartEnzymes™



GENOVIS

INSTRUCTIONS FOR PRODUCT

GingisKHAN® 2000 units (B0-GKH-020)

Digestion of up to 2 mg human IgG1

1 Prepare GingisKHAN®

- Reconstitute GingisKHAN in 200 μ l ddH₂O to a concentration of 10 units/ μ l.
- Reconstitue GingisKHAN Reducing Agent in 50 μ L ddH₂O¹.

**2 Add GingisKHAN®**

- Add 1 unit GingisKHAN / 1 μ g IgG.
- Add 1/10 (v/v) freshly prepared GingisKHAN Reducing Agent.

**3 Digestion**

Incubate for 1-2 h² at 37°C.



PRODUCT DESCRIPTION

GingisKHAN (Gingipain K) is an enzyme that digests human IgG1, yielding intact Fab and Fc fragments. In native conditions, GingisKHAN digests human IgG1 at one single site in the upper hinge (Table 1).

A second digestion site on the Fc may appear if the N-glycans are removed. GingisKHAN is a cysteine protease and requires reducing conditions to be active. Intact Fab and Fc fragments are obtained with GingisKHAN digestion of human IgG1, since mild reducing conditions (i.e. 2 mM cysteine) is sufficient for enzyme activity. Optimal activity is obtained at 37°C and pH 8. The reducing agent is supplied together with the enzyme.

GingisKHAN is purified from *Porphyromonas gingivalis*.

Table 1. *GingisKHAN digestion site.*

| IgG subclass | GingisKHAN digestion site |
|---------------------|----------------------------------|
| Human IgG1 | ..KSCDK / THTCPPCP.. |

Unit Definition

One unit GingisKHAN digests $\geq 95\%$ of 1 μg human IgG1 when incubated in 0.1 M Tris, pH 8.0 at 37°C for 1 h.

Content and Storage

GingisKHAN is supplied lyophilized in 100 mM Tris, 75 mM NaCl, pH 8.0, with no preservatives added. It is supplied with 5 vials of lyophilized GingisKHAN Reducing Agent (10x), yielding 20 mM cysteine upon reconstitution (neutral pH).

GingisKHAN and GingisKHAN Reducing Agent are shipped cold and should be stored at -20°C upon arrival.

After reconstitution, GingisKHAN is stable for 2 months at +4-8°C. Reconstituted GingisKHAN Reducing Agent should be used the same day as it is prepared, it cannot be stored.

GingisKHAN is for R&D use only.

Additional Materials Supplied

- GingisKHAN Reducing Agent (10x), 5 vials.

Additional Materials Required

- Digestion buffer³: 0.1 M Tris, pH 8.0.

Preparations

- Prepare the IgG (human IgG1) in digestion buffer. The final IgG concentration should be 0.5-5 mg/ml.
- GingisKHAN Reducing Agent must be prepared and used the same day, it cannot be stored¹.

Generation of Fab fragments from hIgG1 with GingisKHAN is adversely affected at denaturing conditions i.e. in the presence of chaotropic agents and/or detergents. If the analysis of the digestion efficiency is done with SDS-PAGE, stop the digestion reaction by adding 10 mM iodoacetamide before SDS is added to the SDS-PAGE sample preparation.

Antibody Subunit Generation

1 Prepare GingisKHAN®

- Reconstitute GingisKHAN in 200 μ l ddH₂O to a concentration of 10 units/ μ l.
- Reconstitute GingisKHAN Reducing Agent in 50 μ l ddH₂O and keep on ice¹.
Note! Use the same day as it is prepared, it cannot be stored.

2 Add GingisKHAN®

- Add 1 unit GingisKHAN / 1 μ g IgG.
- Add GingisKHAN Reducing Agent to the reaction mixture. Add 1/10 (v/v) to yield 2 mM cysteine in the reaction.

3 Digestion

- Incubate for 1-2 hours² at 37°C.

Notes

1. Upon reconstitution, the GingisKHAN Reducing Agent may appear cloudy. This does not affect the performance. Make sure to mix it thoroughly before using it in the reaction.
2. Digestion time may need to be optimized for individual antibodies.
3. Other buffers at pH 7-8 can be used, but optimization is required. Sodium chloride concentrations above 75 mM may negatively affect enzymatic activity.

Quality Control

GingisKHAN is tested to meet the specifications and lot-to-lot consistency.

GingisKHAN is tested for absence of microbial contamination with blood agar plates, Sabouraud dextrose agar plates and fluid thioglycollate medium.

Related Products

GingisKHAN® Fab kit

Generation and purification of Fab fragments from hlgG1

FabALACTICA™

Generation of Fab fragments from hlgG1

GingisKHAN®

Limited Use Label License: Research Use Only

All rights reserved. The trademark GingisKHAN® is the property of Genovis AB. GingisKHAN® is duly registered in EU, Canada, USA, Norway, Singapore, China, Australia and Japan.

For research use only. Not intended for any animal or human therapeutic or diagnostic use.

All goods and services are sold subject to Genovis' General Terms and Conditions of Sale.


©2017 Genovis AB

FragIT™ kit

Digestion of IgG and Purification of F(ab')₂ and Fc Fragments

FragIT kit consists of an IgG digestion column, FragIT, and an affinity purification column, CaptureSelect™*. FragIT is a resin with FabRICATOR® enzyme covalently coupled to agarose beads for subunit fragmentation of IgG to generate F(ab')₂ and Fc fragments. After digestion, the fragments can be purified using the CaptureSelect column supplied in the kit.

- Digestion of IgG on a column.
- Purification of F(ab')₂ and Fc fragments.

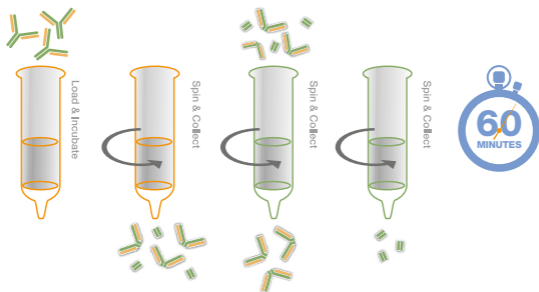


**Made with Thermo Scientific™ CaptureSelect™ resin from Thermo Fisher Scientific Inc. and its subsidiaries. Thermo Scientific and CaptureSelect are trademarks of Thermo Fisher Scientific Inc. and its subsidiaries.*

GingisKHAN[®] Fab kit

Generation and Purification of Fab Fragments from Human IgG1

GingisKHAN Fab Kit consists of GingisKHAN enzyme for upper hinge digestion of human IgG1 and an affinity column for purification of Fab fragments. After digestion, the Fab fragments can easily be purified using the CaptureSelect™ CH1* specific spin column supplied in the kit.

- Digestion of human IgG1 above the hinge.
- Purification of intact Fab fragments.

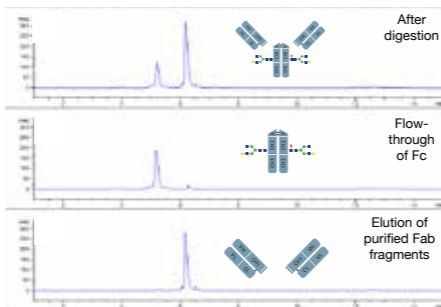

**Made with Thermo Scientific™ CaptureSelect™ resin from Thermo Fisher Scientific Inc. and its subsidiaries. Thermo Scientific and CaptureSelect are trademarks of Thermo Fisher Scientific Inc. and its subsidiaries.*



GENOVIS

US & Canada

Genovis Inc.
245 First Street, Suite 1800
Cambridge, MA 02142
USA

Customer service: 001 (617)-444-8421
Order phone (toll free): 001 (855)-782-0084
Order fax: 001 (858)-524-3006
Email: orders.us@genovis.com

EMEA & Asia

Genovis AB
Box 790
SE-220 07 Lund
Sweden

Customer service: 0046 (0)46 10 12 30
Order phone: 0046 (0)46 10 12 30
Order fax: 0046 (0)46 12 80 20
Email: order@genovis.com