



# OpeRATOR™

Lyophilized

STORE AT

**-20°C**



FOR RESEARCH USE ONLY

## Instructions for Use

OpeRATOR™ Lyophilized 2000 units (G2-OP1-020)  
Process 2 mg O-glycoprotein





## Preparations

### Additional Materials Required

- Reaction buffer: 20mM Tris, pH 6.8.<sup>2</sup>

2. OpeRATOR has optimal activity in pH 5.5-7.5.

## O-glycan-specific Protein Digestion

### Sample Preparation

Prepare the glycoprotein in the reaction buffer to a concentration of 0.1-2 mg/ml.

#### 1. Prepare OpeRATOR and SialEXO

1.1 Reconstitute OpeRATOR and SialEXO in 50  $\mu$ l ddH<sub>2</sub>O each, to a concentration of 40 units/ $\mu$ l.

#### 2. Add SialEXO

2.1 Add 1 unit SialEXO / 1  $\mu$ g glycoprotein.<sup>3</sup>

#### 3. Add OpeRATOR

3.1 Add 1 unit OpeRATOR / 1  $\mu$ g glycoprotein.<sup>3</sup>

#### 4. Enzymatic Reaction

4.1 Incubate for 2 h to overnight (16-18h) at 37°C.

Note: The activity of the OpeRATOR enzyme is significantly reduced if sialic acids are present.<sup>1</sup>

3. A higher enzyme concentration may increase digestion efficiency of individual glycoproteins. This requires optimization.



**USA & Canada**

Genovis Inc.

10919 Technology Place Suite C, San Diego, CA 92127, USA

Phone: 1-855-782-0084 (toll free)

Fax: 1-858-524-3006

**EMEA & Asia**

Genovis AB

Box 4, SE-24421 Kävlinge, Sweden

Phone: +46 46 10 12 30

Fax: +46 46 12 80 20

[support@genovis.com](mailto:support@genovis.com)

[www.genovis.com](http://www.genovis.com)



All rights reserved. Genovis products may be covered by one or more patents, trademarks and copyrights owned or controlled by Genovis AB. For more information about commercial rights, please contact the Genovis team at [licensing@genovis.com](mailto:licensing@genovis.com).

Genovis products are intended for research use only. They are not intended to be used for therapeutic or diagnostic purposes in humans or animals.

All goods and services are sold subject to Genovis' General Terms and Conditions of Sale.

© Genovis AB