



PNGase F

Lyophilized

STORE AT

-20°C



FOR RESEARCH USE ONLY

Instructions for Use

PNGase F Lyophilized 1000 units (G1-PF1-010)

Process 1 mg glycoprotein

PNGase F Lyophilized 5 × 1000 units (G1-PF1-050)

Process 5 × 1 mg glycoprotein

DOWNLOAD INSTRUCTIONS FOR USE



www.genovis.com/ifu-G1-PF1

Lyophilized Enzyme for Hydrolysis of N-glycans on Glycoproteins

PNGaseF (Peptide N-glycosidase F) is a glycoamidase hydrolyzing the amide bond between the polypeptide asparagine and the innermost GlcNAc of all mammalian asparagine-linked complex, hybrid, or high-mannose oligosaccharides. The activity of PNGaseF on some glycoproteins can be slow or inhibited due to steric hindrance - longer incubation times or denaturation of the glycoprotein may in these cases be required.

PNGaseF is derived from *Elizabethkingia meningoseptica* and expressed in *E. coli*. The enzyme contains a His-tag and has a molecular weight of 36 kDa.

UNIT DEFINITION

1 unit PNGaseF Lyophilized removes N-glycans from $\geq 95\%$ of 1 μg glycoprotein (etanercept) when incubated in PBS (10 mM sodium phosphate, 150 mM NaCl) pH 7.4 at 37°C for 1 hour.

CONTENT AND STORAGE

PNGase F Lyophilized is supplied lyophilized in 50 mM HEPES buffer pH 7.5, with no preservatives added.

PNGase F Lyophilized is shipped cold and should be stored at -20°C upon arrival. After reconstitution, PNGaseF Lyophilized is stable for at least 1 month at +4-8°C.

PNGase F Lyophilized is for R&D use only.

QUALITY CONTROL

PNGaseF Lyophilized is tested to meet the specifications and lot-to-lot consistency.

PNGaseF Lyophilized is tested for absence of microbial contamination with blood agar plates, Sabouraud dextrose agar plates and fluid thioglycollate medium.

YOU MIGHT ALSO BE INTERESTED IN

PNGase F Immobilized

Immobilized enzyme for hydrolysis of N-glycans in spin columns

OglyZOR™

Hydrolysis of core 1 O-glycans

GlycINATOR™

Hydrolysis of all Fc N-glycans

OmniGLYZOR®

Hydrolysis of N- and mucin-type O-glycans

Preparations

Additional Materials Required

- Reaction buffer: PBS pH 7.4.¹

1. Other common buffers in pH 6.0-9.0 over a wide range of ionic strengths (0-1000mM NaCl) may also be used. Optimization may be required if buffers other than the recommended are used.

Hydrolysis of N-glycans on Glycoproteins

Sample Preparation

Prepare the glycoprotein in the reaction buffer. The final glycoprotein concentration in the reaction should be 0.5-5 mg/ml.

1. Prepare PNGase F

1.1 Reconstitute PNGase F in 50 μ l ddH₂O to a concentration of 20 units/ μ l.

2. Add PNGase F

2.1 Add 1 unit PNGase F / 1 μ g glycoprotein.²

3. Enzymatic Reaction

3.1 Incubate for 1 h to overnight at 37°C.³

Note: Optimization of the enzyme concentration, incubation time and incubation temperature may be needed depending on the substrate. Some glycoproteins require denaturing and/or reducing reaction conditions to reach full deglycosylation.

2. A higher enzyme concentration may increase digestion efficiency of individual glycoproteins. This requires optimization.
3. Incubation temperature can be increased to up to 50°C for a faster deglycosylation. A surfactant may be required to prevent precipitation of the substrate protein.

USA & Canada

Genovis Inc.

10919 Technology Place Suite C, San Diego, CA 92127, USA

Phone: 1-855-782-0084 (toll free)

Fax: 1-858-524-3006

EMEA & Asia

Genovis AB

Box 4, SE-24421 Kävlinge, Sweden

Phone: +46 46 10 12 30

Fax: +46 46 12 80 20

support@genovis.com

www.genovis.com



All rights reserved. Genovis products may be covered by one or more patents, trademarks and copyrights owned or controlled by Genovis AB. For more information about commercial rights, please contact the Genovis team at licensing@genovis.com.

Genovis products are intended for research use only. They are not intended to be used for therapeutic or diagnostic purposes in humans or animals.

All goods and services are sold subject to Genovis' General Terms and Conditions of Sale.

© Genovis AB