



FabRICATOR®Z

Lyophilized

STORE AT

-20°C



FOR RESEARCH USE ONLY

Instructions for Use

FabRICATOR®Z Lyophilized 2000 units (A0-FRZ-020)
Process 2 mg mouse IgG2a and IgG3

DOWNLOAD INSTRUCTIONS FOR USE



www.genovis.com/ifu-A0-FRZ

Lyophilized Enzyme for Below Hinge Digestion of Mouse IgG2a and IgG3

FabRICATOR Z (IdeZ) is a cysteine protease that digests mouse IgG2a and IgG3 at a single amino acid site below the hinge, generating homogenous F(ab')₂ and Fc fragments. There is no risk of overdigestion if the incubation time is prolonged. Since FabRICATOR Z digests IgG under physiological reaction conditions, the immunoreactivity is preserved. Most mouse IgG2a and IgG3 are digested by FabRICATOR Z within 2 hours, although the enzyme:antibody ratio and/or digestion time may need to be optimized for certain antibodies. FabRICATOR Z is also active on IgG from monkey, rabbit and sheep. For digestion of human IgG, we recommend using FabRICATOR (A0-FR1-020).

FabRICATOR Z Lyophilized is derived from *Streptococcus equi subsp. zooepidemicus* and expressed in *E. coli*. The enzyme contains a His-tag and the molecular weight is 36 kDa.

UNIT DEFINITION

One unit FabRICATOR Z Lyophilized digests $\geq 95\%$ of 1 μg mouse IgG2a when incubated in 10 mM sodium phosphate, 10 mM NaCl, pH 6.5 at 37°C for 2 hours.

CONTENT AND STORAGE

FabRICATOR Z Lyophilized is supplied lyophilized in 50 mM Tris, 150 mM NaCl, pH 7.6, with no preservatives added.

FabRICATOR Z Lyophilized is shipped at ambient temperature and should be stored at -20°C upon arrival.

After reconstitution, FabRICATOR Z Lyophilized is stable for at least 1 month at +4-8°C.

FabRICATOR Z Lyophilized is for R&D use only.

QUALITY CONTROL

FabRICATOR Z Lyophilized is tested to meet the specifications and lot-to-lot consistency.

FabRICATOR Z Lyophilized is tested for absence of microbial contamination with blood agar plates, Sabouraud dextrose agar plates and fluid thioglycollate medium.

YOU MIGHT ALSO BE INTERESTED IN

FabRICATOR® Z Immobilized

Immobilized enzyme for below hinge digestion of mouse IgG2a and IgG3 in spin columns

FabRICATOR® Z Fab2 Kit

Immobilized enzyme and affinity resin for below hinge digestion of mouse IgG2a and IgG3 and purification of fragments

FabRICATOR® Lyophilized

Lyophilized enzyme for below hinge digestion of IgG

FabALACTICA® Lyophilized

Lyophilized enzyme for above hinge digestion of human IgG1

Preparations

Additional Materials Required

- Reaction buffer: 10 mM sodium phosphate, 10 mM NaCl, pH 6.5.¹

1. The NaCl concentration and pH can be increased but this may require increased enzyme concentration and/or a prolonged incubation time.

Below Hinge Digestion of Mouse IgG2a and IgG3

Sample Preparation

Prepare the IgG in the reaction buffer. The final IgG concentration in the reaction should be 0.5-10mg/ml.

1. Prepare FabRICATOR Z

1.1 Reconstitute FabRICATOR Z in 50 μ l ddH₂O to a concentration of 40 units/ μ l.

2. Add FabRICATOR Z

2.1 Add 1 unit FabRICATOR Z / 1 μ g IgG.²

3. Enzymatic Reaction

3.1 Incubate for 2 h³ at 37°C.

2. The enzyme:antibody ratio may need to be increased for individual antibodies.
3. The incubation time may need to be prolonged for individual antibodies.

USA & Canada

Genovis Inc.

245 First Street, Suite 1800, Cambridge, MA 02142, USA

Phone: 1-855-782-0084 (toll free)

Fax: 1-858-524-3006

EMEA & Asia

Genovis AB

Box 4, SE-24421 Kävlinge, Sweden

Phone: +46 46 10 12 30

Fax: +46 46 12 80 20

support@genovis.com

www.genovis.com

All rights reserved. Genovis products may be covered by one or more patents, trademarks and copyrights owned or controlled by Genovis AB. For more information about commercial rights, please contact the Genovis team at licensing@genovis.com.

Genovis products are intended for research use only. They are not intended to be used for therapeutic or diagnostic purposes in humans or animals.

All goods and services are sold subject to Genovis' General Terms and Conditions of Sale.

© Genovis AB